

ClearH₂O[®]

AquaValve[®]

Drinking Valve for AquaPak[®]

AquaValve[®] is the drinking valve developed for exclusive use with AquaPak[®] water pouches.

Delivering Solutions

- Easy activation for weanling transition
- Durable design for extended use
- Robust, chew-resistant construction
- Convenient bulk packaging
- Fits pouch-compatible and disposable cage systems

Delivering Quality

- Ethylene oxide treated
- FDA-compliant for food and water contact surfaces
- Made from BPA-free POM material
- Certificate of analysis by manufactured lot




Ordering Information

AquaValve[®]
90-07
100 count pack
18 packs/case



ClearH₂O[®]

www.clearh2o.com

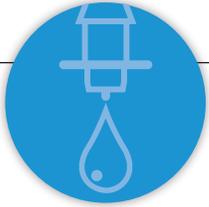

AquaValve[®]

Using AquaValve[®]



1 Remove the cage lid and align the AquaValve[®] for insertion into the lid water access hole in the water bottle receptacle.



2 Support the bottom of the lid with two fingers on either side of the water access hole as you push the activator end of the valve through the hole.



3 With a relaxed grip, orient AquaPak[®] with the smooth side approaching the piercing cone of the AquaValve[®]. Keep two fingers on the bottom of the lid to support the valve.



4 With a firm grip to pressurize the bottom of the AquaPak[®], thrust the pouch forward piercing it onto the AquaValve[®] collar. Check the AquaValve[®] by gently touching the activator to observe water flow.

ClearH₂O[®]

Company Information

ClearH₂O[®] is committed to advancing animal research and agriculture with more effective and ethical hydration, nutrition, and medication delivery solutions utilizing our proprietary gel technology. Made in our FDA approved facility, ClearH₂O[®] products are formulated with purified ingredients and highly palatable to optimize consumption. The ready to use, convenient product format with long-dated, room temperature shelf stability helps animal researchers and producers improve the health and welfare of their animals while optimizing research and productivity outcomes.

www.clearh2o.com

85 Bradley Drive Westbrook, ME 04092 1-888-HYDROGL (1-888-493-7645) Fax: 1-207-774-3449